

CELLMATE® II

A Lifetime of Power



The World's Most Advanced Serological Pipettor

CellMate II is a uniquely designed lightweight, cordless pipettor for use with graduated and volumetric glass and plastic serological pipettes. Ergonomically designed, CellMate II offers simple efficient pipetting performance and maximum pipetting comfort.

Featuring a heavy-duty rechargeable lithium ion battery that's guaranteed for life, this pipettor can operate approximately 3 to 4 times longer between charges than typical units using nickel cadmium batteries.

Available in a choice of five colors, the CellMate II can be wall- or table-mounted. A 1mL pipette support is included to eliminate wobble when using 1mL pipettes.

The CellMate II is simply the best cordless pipettor you can buy!

FEATURES

- Lifetime guaranteed lithium ion battery
- Lightweight ergonomic design
- Separate aspirate and dispense speed controls
- Fully autoclavable nosecone housing
- Large backlit LED display
- Customizable ID area, 5 color options
- 1mL pipette support

BENEFITS

- Offers longer usable power per charge, more charges and no battery memory
- Allows longer, fatigue-free pipetting
- Affords greater control over procedures
- Facilitates disinfection and decontamination
- Provides visual inspection of remaining battery charge and speed settings
- Permits easy recognition of ownership and/or task designation
- Creates additional support when using 1mL pipettes


fluid thinking.
www.matrixtechcorp.com

Cordless Pipettor With Cutting Edge Battery Technology

Special circuitry on the CellMate II makes it impossible to overcharge the battery and allows the pipettor to be charged or partially charged at any convenient time. A battery icon on the large backlit LCD display indicates the level of charge, and the display blinks when the battery charge is getting low.



Buttons on the rear of the CellMate II unit allow you to set aspiration (upper button) and dispense (lower button) speeds. Eight speeds are available for each function – simply push “+” for faster and “-” for slower. Your selection will be displayed on the lower portion of the pipettor’s LCD. A zero speed selection on dispense enables gravity dispensing.



The entire CellMate II can be cleaned with most common laboratory disinfectants. The fully autoclavable nosecone housing contains a silicone rubber pipet holder, a 0.45 micron hydrophobic filter assembly and a special 1mL pipette-stabilizing attachment.



A customizable ID area on the CellMate II unit allows individual user names to be displayed on the body of the pipettor for easy identification. As an added means of differentiation, the CellMate II is offered in five color options.



The CellMate II kit includes both a table-mounted stand and a wall-mounted holder.



CellMate II

Item No.	Description	Unit
9501	Cellmate II (White)	
9511	Cellmate II (Clear)	
9521	Cellmate II (Blue)	
9531	Cellmate II (Red)	
9541	Cellmate II (Green)	
S▶ 9056	Hydrophobic Filters 0.8µm	24/case
S▶ 9057	Hydrophobic Filters 0.45µm	24/case
9064	Nosepiece (holder, silicone gripper, filter)	
9065	Silicone pipet gripper	4/case
9066	Wall Mount	
9067	Stand	
9068	Power Supply	
9070	1mL Pipet Support	

S▶ =Sterile

ORDERING GUIDE

HANDHELD PIPETTING

AUTOMATED PIPETTING

CONSUMABLES

Corporate Headquarters
22 Friars Drive
Hudson, NH 03051 USA
Tel: 603-595-0505
800-345-0206
e-mail: matrix@matrixtechcorp.com

European Headquarters
Lower Meadow Road
Brooke Park
Handforth, Wilmslow
Cheshire, SK9 3LP, UK
Tel: +44 (0) 161 486 2110
e-mail: matrix@matrixtechcorp.eu.com

Asia-Pacific Headquarters
Akemi-14-11-705
Urayasu-City
Chiba-Pref
279-0014 Japan
Tel: 81 (0) 47 355 6491
e-mail: asia@matrixtechcorp.com

MATRIX
fluid thinking.

www.matrixtechcorp.com