
Biohit Midi Pro Operator Instructions

Pipette Controller

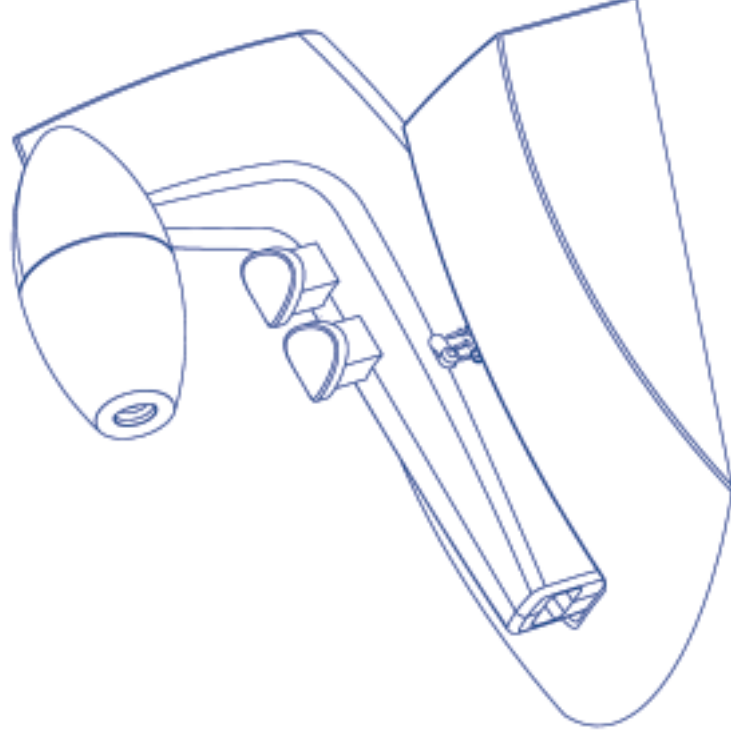BIOHIT

Operator Instructions

2 – 8

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9 – 15

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
16 – 22

Manual Del Operador

23 – 29

Notes

30 – 31

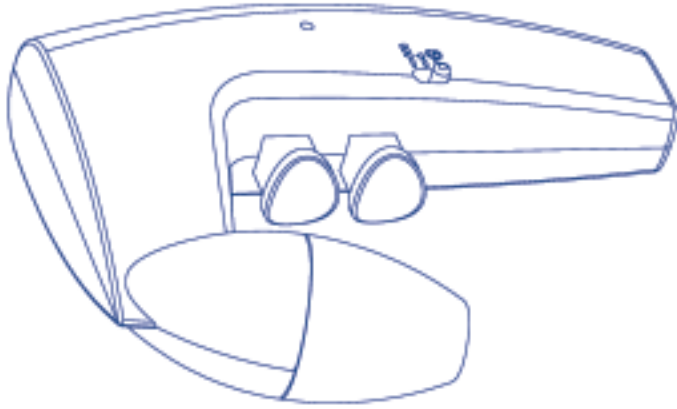<p>Biohit Plc. Headquarters Laippatie 1, 00880 Helsinki, Finland Tel: +358-9-773 861 Fax: +358-9-773 86200 E-mail: info@biohit.com www.biohit.com</p>	<p>CHINA Finland Biohit Co., Ltd. Tel: +86-21-6248 5589 Fax: +86-21-6248 7786 E-mail: info.china@biohit.com</p>	<p>FRANCE Biohit S.A. Tel: +33-1-3088 4130 Fax: +33-1-3088 4102 E-mail: commercial.france@biohit.com</p>
<p>GERMANY Biohit Deutschland GmbH Tel: +49-6003-82 820 Fax: +49-6003-82 8222 E-mail: info@biohit.de</p>	<p>JAPAN Biohit Japan Co., Ltd. Tel: +81-3-5822 0021 Fax: +81-3-5822 0022 E-mail: sales@biohit.co.jp</p>	<p>RUSSIA Biohit OOO. Tel: +7-812-327 5327 Fax: +7-812-327 5323 E-mail: main@biohit.ru</p>
<p>U.K. Biohit Ltd. Tel: +44-1803-315 900 Fax: +44-1803-315 530 E-mail: info@biohit.co.uk</p>	<p>U.S.A. Biohit Inc. Tel: +1-732-922-4900 Fax: +1-732-922-0557 E-mail: pipet@biohit.com</p>	<p>CE BIOHIT</p>

Introduction

The Biohit Midi Pro has been carefully designed and manufactured as a lightweight, cordless pipetting aid for use with all plastic or glass pipettes from 1–100ml.

This new Midi Pro is constructed in recyclable materials and operates on two rechargeable, long life, environmentally friendly, 1.2V nickel metal hydride batteries. Biohit Midi Pros have uniquely formed finger triggers to minimize any high pressure points. These triggers also control intake and dispense speeds.


A battery warning light (4) is triggered whenever the charge in the batteries is low and they should then be recharged. Pipetting with Biohit Midi Pros can also take place while the units are connected to their chargers (5).

Midi Pro have a conveniently located slide switch (3) which allows the user to select between three modes of operation:

Slide switch position 'High' (top)	Suction and dispense at full speed (Blow-out)
Slide switch position 'Low' (middle)	Suction and dispense at reduced speed (Blow-out)
Slide switch position 'G' (bottom)	Suction at reduced speed, dispense by gravity (To Deliver)

Technical Data

Batteries	2 x 1.2V nickel metal hydride
Battery service Life	500–1000 cycles from empty
Charger	75mA/h for 14 hours plug-in; UK, European, USA, Japanese or Australian versions
Overall weight	Installation (over voltage) category II
Filter (8)	Double insulation indicated on charger by symbol 
Case material	180g without stand
Nosecone material	Hydrophobic 2.5mm dia 0.45µm (fitted) plus two additional spare filters
Pipette holder	1x 0.45µm
Pipette types	1x 0.20µm ASA (10) ASA Silicone (9) Plastic or glass 1–100ml

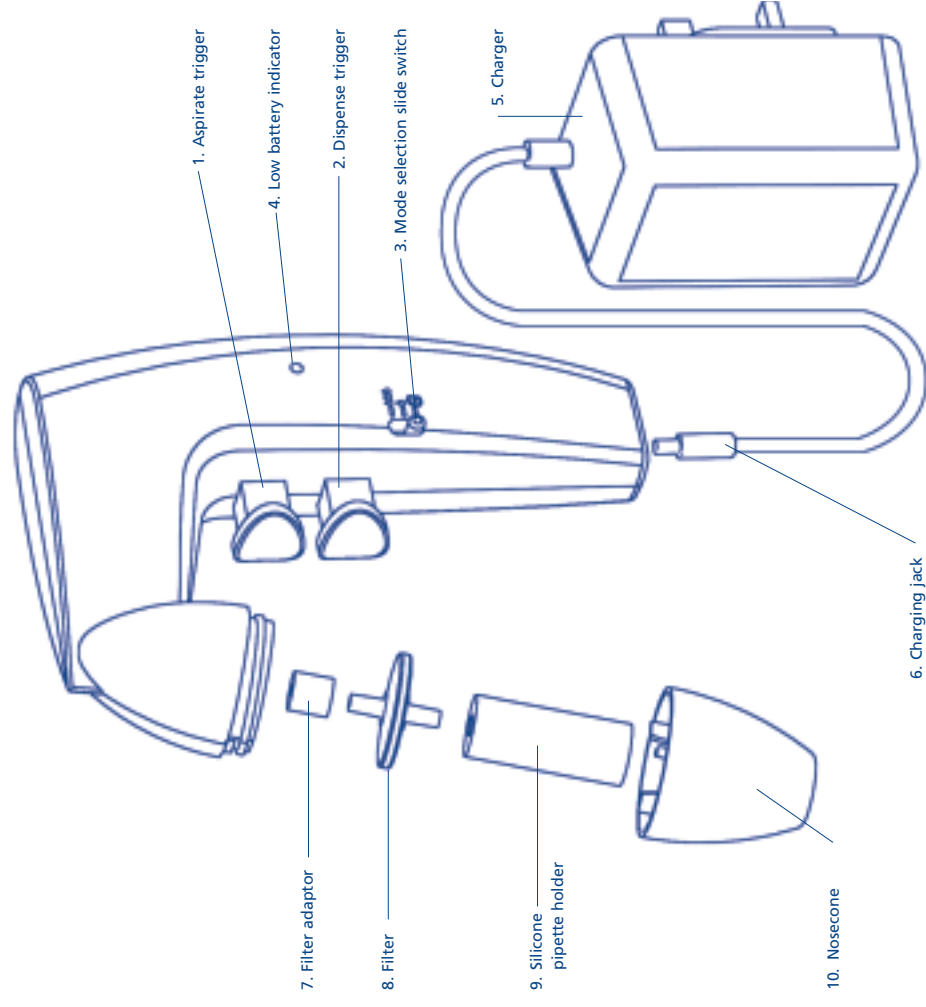Contents

The Midi Pro kit comprises

- (a) Midi Pro pipette controller
- (b) Charger; either
 - 220V-240V 50/60Hz 3-pin (UK),
 - 100V-120V 50/60Hz 2-pin (USA),
 - 220V-240V 50/60Hz 2-pin (Australia)
- (c) Two spare hydrophobic filters, 1 x 0.45µm and 1 x 0.20µm
- (d) Bench stand / wall bracket
- (e) Instruction manual

Safety Instructions

The Midi Pro has been designed with safety in mind. They are lightweight and balance safely in the hand. The triggers have been carefully designed to eliminate pressure points and control of intake and delivery is therefore made easier.

It is recommended when using deleterious liquids that the user wears protective clothing including goggles and gloves. If in doubt ask your safety officer. Please consult and establish appropriate health and safety practices applicable to the liquid being dispensed within your working environment.

WARNING!

- Only use this instrument within the recommended limits of its chemical resistance.
- Never use with fuming liquids which are incompatible with ASA, ABS, silicone or polypropylene.
- Never use with highly flammable liquids.
- Observe the reagent manufacturer's safety information.
- Never use the instrument in an atmosphere with danger of explosion.
- When not in use store on bench stand or wall bracket.
- Never use force when in use.
- Should the instrument fail to work to specification stop pipetting and check instructions headed Trouble Shooting (page 7).
- Only recharge with manufacturer's original charger.
- All repairs must be carried out by an authorized dealer.
- Only use the manufacturer's accessories and spare parts.
- Charger must only be used indoors and protected against moisture.

Recharging the Midi Pro

Before starting to operate, carefully check mains voltage and recharge batteries.

To charge the Midi Pro


- 1 Recharging temperature +10°C to +35°C.
 - 2 Plug the charger jack (6) into the base of the Midi Pro.
 - 3 Plug the charger (5) into a suitable wall socket. The red indicator on the charger will light up.
 - 4 Place Midi Pro onto bench stand.
 - 5 Allow unit to charge for 14 hours from empty to fully charged.
 - 6 Once the Midi Pro is fully charged, the charger automatically switches to a trickle charge operation and the red indicator will turn off. If the charge to the Midi Pro is interrupted and then re-started, the charger starts its full 14 hour charging cycle again.
- This will not damage the battery.

Only use the charger indoors and protect against moisture.


Fitting the Midi Pro wall bracket

The Midi Pro bench stand is supplied ready to use as a wall bracket. When used as a bench stand it is recommended that the wall fitting is removed.

Bench stand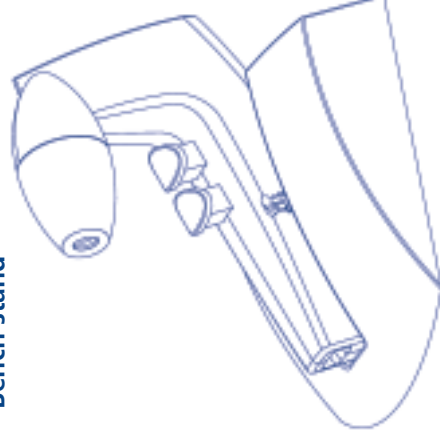

Wall bracket

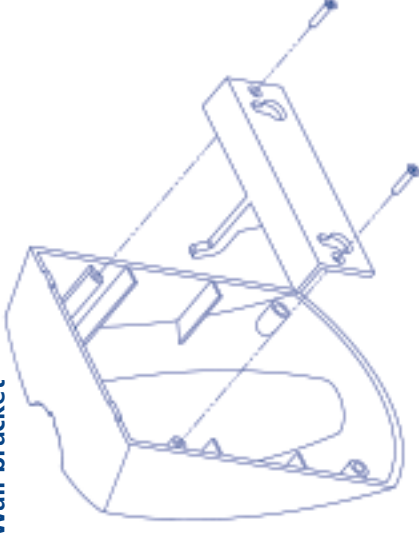
Cleaning

Before cleaning disconnect from the mains charger. Do not clean the mains charger.

The Midi Pro can be cleaned with most common laboratory disinfectants. The silicone pipette holder (9) can be autoclaved.

To remove the nosecone and pipette holder turn the nosecone anti-clockwise. When the nosecone is removed the silicone pipette holder can also be disengaged from the filter by gently turning.

The hydrophobic filter (8) will require occasional replacement when wetted by the liquid and whenever suction capacity is slowing.

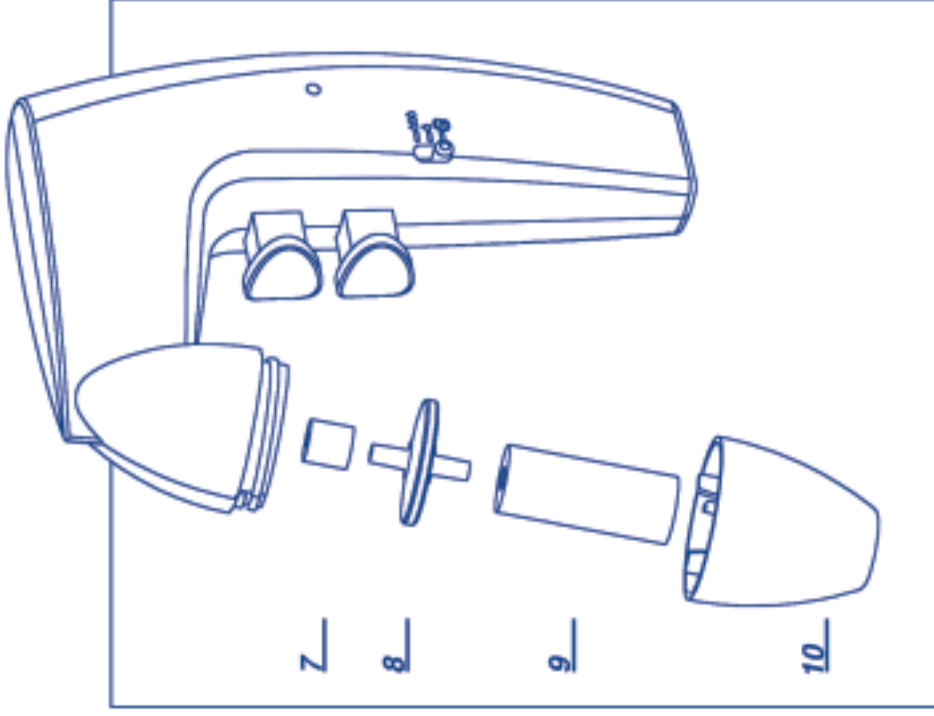Pipetting

To aspirate

- 1 Carefully hold the pipette and insert it securely into the polypropylene nose cone and silicone holder.
- 2 Select mode of operation—High, Low, for blowout pipettes. 'G' for 'to deliver'.
- 3 Depress upper trigger (1) (The harder the pipetting triggers are pressed the faster the aspiration.)

To dispense

- 1 In BLOW-OUT mode, (mode switch on high or low) depress the lower trigger (2) (The harder the pipetting triggers are pressed, the faster the dispense).
- 2 If 'G' gravity mode is chosen, the upper aspirate trigger automatically operates at low speed while the lower trigger opens the pipette to atmosphere, allowing liquid to run out.

WARNING!

Please ensure that no liquid enters the instrument. The Midi Pro should never be used without the filter (8). The filter also prevents liquid from entering the instrument. Where liquid has entered please clean and replace filter.

Trouble Shooting

The Midi Pro is virtually service-free. If the recommendations below do not resolve the problem, please contact your authorized dealer.

TROUBLE	POSSIBLE CAUSE	REMEDY
Suction capacity slow	Filter blocked, dirty or damaged	Clean or replace membrane filter
	Batteries low	Recharge
Pipette drips	Pipette not securely positioned in silicone pipette holder	Re-attach
	Pipette holder or pipette damaged	Replace

Note

When the meniscus level has been reached liquid should not drip out of the pipette. If the pipette does drip, check that the adaptor assembly and filter are correctly assembled and the pipette is properly located into the silicone pipette holder.

Return for Repair

In the unlikely event of repair, or where the damage to the unit necessitates return to your distributor or the manufacturer, please decontaminate the unit and complete the required decontamination certificate where appropriate. Please also give details of the fault.

Voltages

UK	220V-240V	50/60Hz	3-pin
Europe	220V-240V	50/60Hz	2-pin
USA	100V-120V	50/60Hz	2-pin
Japan	100V-120V	50/60Hz	2-pin
Australia	220V-240V	50/60Hz	2-pin

Ordering Data

Midi Pro

Midi Pro complete with charger, bench stand/ wall bracket with two spare membrane filters.

	UK	Europe	USA	Japan	Aust./NZ
Pro	Cat. No.	Cat. No.	Cat. No.	Cat. No.	Cat. No.
	710913	710911/EUR	710912/US	710914/JPN	710915/AUS/NZ

Box of 5 spare 0.45µm filters721962
Box of 5 spare 0.20µm filters721965
0.45µm sterile filter (1 off)721966
Silicone Pipette holder721963
ASA nosecone (white)721967

WEEE

 In compliance with European Directive, WEEE (2002/96/EC) on waste and reducing of hazardous substances in electrical and electronic equipments, this device must not be recycled as unsorted municipal waste. Instead this device must be collected separately in accordance the local recycling regulations. The solid bar used in conjunction with the crossed-out wheeled bin indicates that the product was put on the European market after 13 August 2005.

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