

The Ovation M pipette is a variable volume, air displacement, 2 stroke pipette intended to aspirate and dispense precise fluid volumes.



****The Ovation M
is autoclavable****

Using your Ovation pipette



Ergonomic Handling


1. "Shake hands" position to pick up. Allow body to fill your palm.
2. Rotate adjustable hook to rest comfortably on your forefinger.
3. Keep hand posture relaxed.



To Set Volume

Rotate the plunger to display the desired volume (clockwise to decrease, counter clockwise to increase).

Do not rotate beyond the pipette's volume limits!

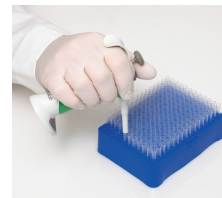

Installing the "Click" Tip

1. Insert nozzle into tip.
2. Lightly press down until a "click" is heard. This indicates the tip is securely sealed onto the nozzle.

Taking Care of Your Ovation Pipette

The Ovation pipette requires minimal routine maintenance. Always store it in its "standing" position when not in use. Clean outer surfaces as needed with a soft cloth dampened with warm water. To decontaminate outer surfaces, wipe with 70% ethanol or isopropanol solution or use a 10% bleach solution followed by water. Remove ejector sleeve and wipe outside of nozzle and inside of ejector sleeve with 70% alcohol to remove any accumulated deposits. **ONLY WIPE THE LCD SCREEN DISPLAY WITH WATER.**

- Only guaranteed to work with VistaLab Tips.
- 2 Year Limited Warranty.
- VistaLab Technologies offers comprehensive calibration and repair programs recommended every 6-12 months.
See www.vistalab.com/services for more information or to place an order.
- For more information on troubleshooting and maintenance visit www.vistalab.com/support.

For additional assistance with warranty repairs or other technical assistance, contact techservice@vistalab.com.

VistaLab Technologies, Inc.
2 Geneva Road
Brewster, NY 10509 USA
(888) 652-6520 or (914) 244-6226

VistaLab™
TECHNOLOGIES