Operator Instructions

1 – 8

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9-16

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17-24

Manual Del Operador

25-32

Jencons Scientific Ltd

Cherrycourt Way Ind. Est.
Stanbridge Rd, Leighton Buzzard
LU7 4UA, ENGLAND

Phone: 01525 372010
Fax: 01525 379547
Int. Phone: +44 1525 372010
Int. Fax: +44 1525 851461

Email: uksales@jencons.co.uk
Int. Email: export@jencons.co.uk
Website: www.jencons.co.uk
Online sales: www.ecomcat.co.uk

Jencons Scientific Inc (USA)

800 Bursca Drive, Suite 801
Bridgeville, PA 15017

Phone: 412 257-8861
Toll Free: 800 846-9959
Fax: 412 257-8809

Email: info@jencons.com
Website: www.jenconsusa.com

ISO 9001:2000

JENCONS


POWERPETTE , POWERPETTE
Plus
& POWERPETTE
turbo

Operator Instructions

Powerpette

Powerpette Plus

Powerpette Turbo

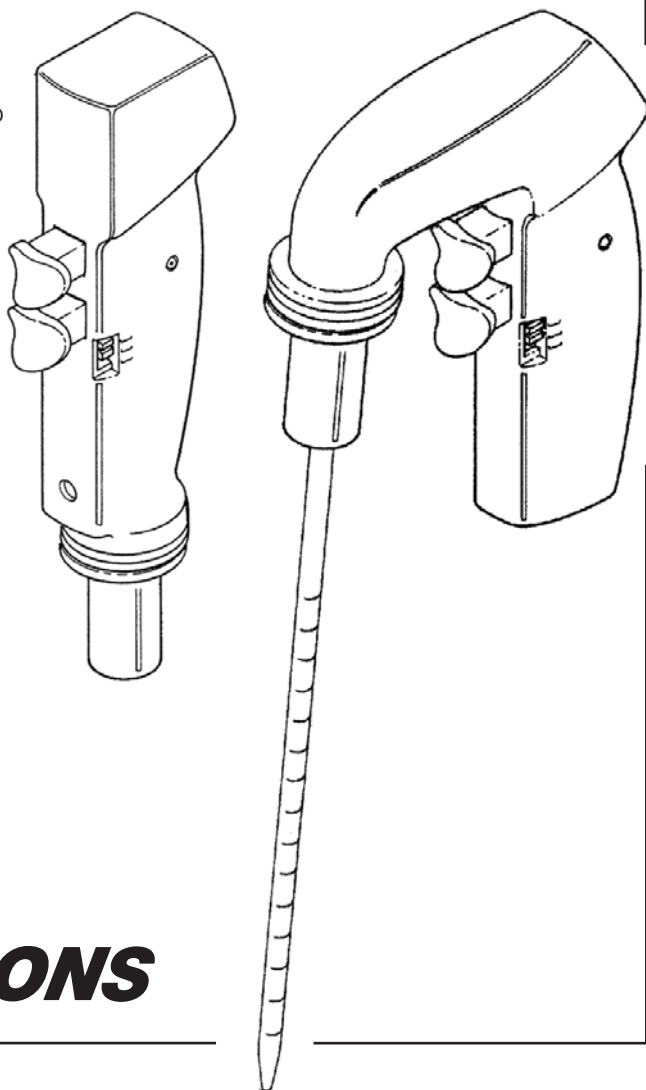JENCONS

Introduction

The Jencons **POWERPETTE**, **POWERPETTE Plus** and **POWERPETTE turbo** have been carefully designed and manufactured as lightweight, cordless pipetting aids for use with all plastic or glass pipettes from 1–100ml.

These new **POWERPETTES** are constructed in recyclable materials and operate on two rechargeable, long life, environmentally friendly, 1.2V nickel metal hydride batteries.

POWERPETTES have uniquely formed finger triggers to minimize any high pressure points. These triggers also control intake and dispense speeds.

A battery warning light (4) is triggered whenever the charge in the batteries is low and they should then be recharged. Pipetting with **POWERPETTES** can also take place while the units are connected to their chargers (5).


POWERPETTES have a conveniently located slide switch (3) which allows the user to select between three modes of operation:

Slide switch position 'High' (top)	Suction and dispense at full speed (Blow-out)
Slide switch position 'Low' (middle)	Suction and dispense at reduced speed (Blow-out)
Slide switch position 'G' (bottom)	Suction at reduced speed, dispense by gravity (To Deliver)

Technical Data

Batteries	2 x 1.2V nickel metal hydride
Battery service Life	500–1000 cycles from empty
Charger	75mA/h for 14 hours plug-in; UK, European, USA, Indian, Japanese or Australian versions Installation (over voltage) category II Double insulation indicated on charger by symbol 
Overall weight	180 g without stand
Filter (8)	Hydrophobic 25 mm dia 0.45 µm (fitted) plus two additional spare filters 1x 0.45 µm 1x 0.20 µm
Manifold material	ABS
Case material	ASA
Nosecone material	Polypropylene (10)
Pipette holder	Silicone (9)
Pipette types	Plastic or glass 1–100 ml

All reference numbers below are applicable to both
POWERPETTE, **POWERPETTE Plus** and **POWERPETTE turbo**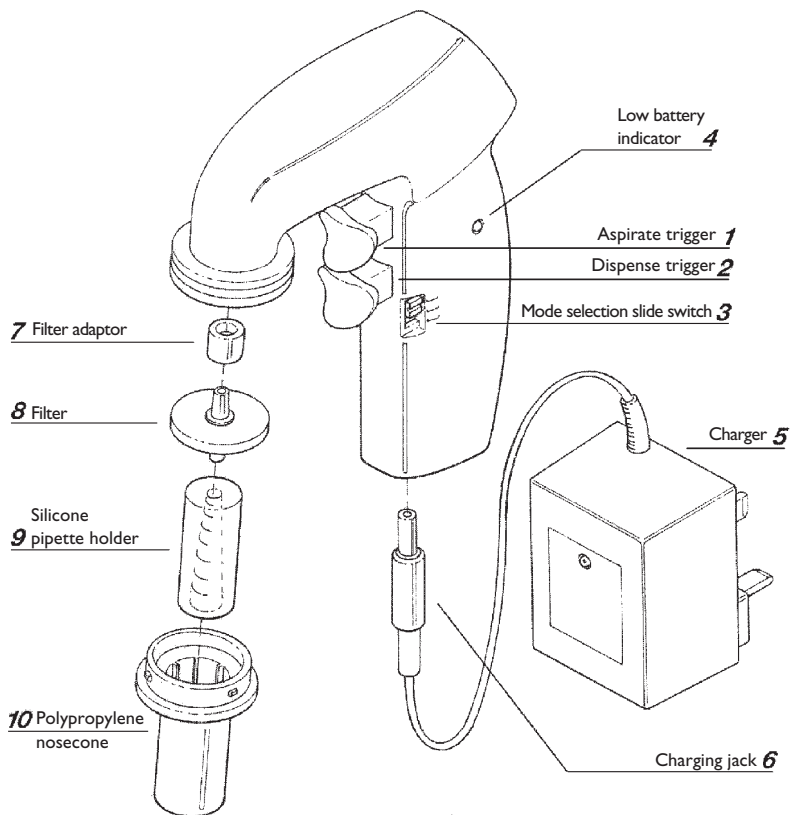

CE **POWERPETTE**, **POWERPETTE Plus** and **POWERPETTE turbo**
are CE marked to Low Voltage Directive 73/23/EEC and
Electromagnetic Compatibility Directive 89/336/EEC

The **POWERPETTE** / **POWERPETTE Plus** / **POWERPETTE turbo** kit comprises

- (a) **POWERPETTE** pipette controller
- (b) Charger; either
220V-240V 50/60Hz 3-pin (UK), 220V-240V 50/60Hz 2-pin (Europe),
100V-120V 50/60Hz 2-pin (USA), 100V-120V 50/60Hz 2-pin (Japan),
220V-240V 50/60Hz 3-pin (India) or 220V-240V 50/60Hz 2-pin (Australia)
- (c) Two spare hydrophobic filters, 1x 0.45 µm and 1x 0.20 µm
- (d) Bench stand (**POWERPETTE** also has a wall bracket)
- (e) Instruction manual

Safety Instructions

The **POWERPETTE**, **POWERPETTE Plus** and **POWERPETTE turbo** have been designed with safety in mind. They are lightweight and balance safely in the hand. The triggers have been carefully designed to eliminate pressure points and control of intake and delivery is therefore made easier.

It is recommended when using deleterious liquids that the user wears protective clothing including goggles and gloves. **If in doubt ask your safety officer.** Please consult and establish appropriate health and safety practices applicable to the liquid being dispensed within your working environment.

WARNING!

- Only use this instrument within the recommended limits of its chemical resistance.
- Never use with fuming liquids which are incompatible with polypropylene or silicone.
- Never use with highly flammable liquids.
- Observe the reagent manufacturer's safety information.
- Never use the instrument in an atmosphere with danger of explosion.
- When not in use store on bench stand or wall bracket.
- Never use force when in use.
- Should the instrument fail to work to specification stop pipetting and check instructions headed Trouble Shooting (page 7).
- Only recharge with manufacturer's original charger.
- All repairs must be carried out by an authorized dealer.
- Only use the manufacturer's accessories and spare parts.
- Charger must only be used indoors and protect against moisture.

Recharging the *POWERPETTE* / *POWERPETTE* Plus / *POWERPETTE* turbo

Before starting to operate, carefully check mains voltage and recharge batteries.

To charge the *POWERPETTE* / *POWERPETTE* Plus / *POWERPETTE* turbo

- 1 Recharging temperature +10°C to +35°C.
- 2 Plug the charger jack (6) into the base of the *POWERPETTE*.
- 3 Plug the charger (5) into a suitable wall socket. The red indicator on the charger will light up.
- 4 Place *POWERPETTE* onto bench stand.
- 5 Allow unit to charge for 14 hours from empty to fully charged.
- 6 Once the *POWERPETTE* is fully charged, the charger automatically switches to a trickle charge operation and the red indicator will turn off. If the charge to the *POWERPETTE* is interrupted and then re-started, the charger starts its full 14 hour charging cycle again.
This will not damage the battery.



Only use the charger indoors and protect against moisture.

Fitting the *POWERPETTE* wall bracket

The *POWERPETTE* wall bracket can be attached by removing the adhesive protective film from the double sided tape and pressing firmly to a grease free smooth surface. Alternatively, it can be attached using the two stainless steel screws provided. Care should be taken in positioning the screws.

Fitting the *POWERPETTE* Plus / *POWERPETTE* turbo into the bench stand

The *POWERPETTE* Plus / *POWERPETTE* turbo can be fitted in the bench stand in one of the two positions shown. It can be charged in either position.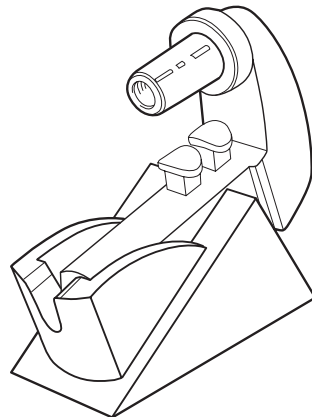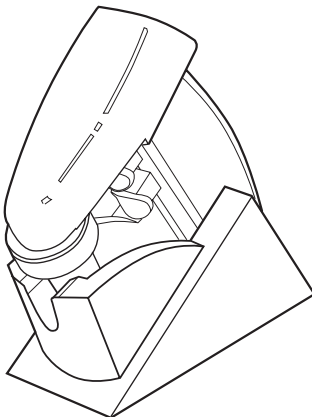

Cleaning

Before cleaning disconnect from the mains charger. Do not clean the mains charger.

The **POWERPETTE** / **POWERPETTE Plus** / **POWERPETTE turbo** can be cleaned with most common laboratory disinfectants. The polypropylene nosecone (10) and silicone pipette holder (9) can be autoclaved.

To remove the nosecone and pipette holder turn the nosecone anti-clockwise. When the nosecone is removed the silicone pipette holder can also be disengaged from the filter by gently turning.


The hydrophobic filter (8) will require occasional replacement when wetted by the liquid and whenever suction capacity is slowing.

Replacing the Filter

The hydrophobic filter (8) will require occasional replacement, particularly when suction capacity is slowing.

Note

Only use genuine manufacturer's recommended filters and accessories.

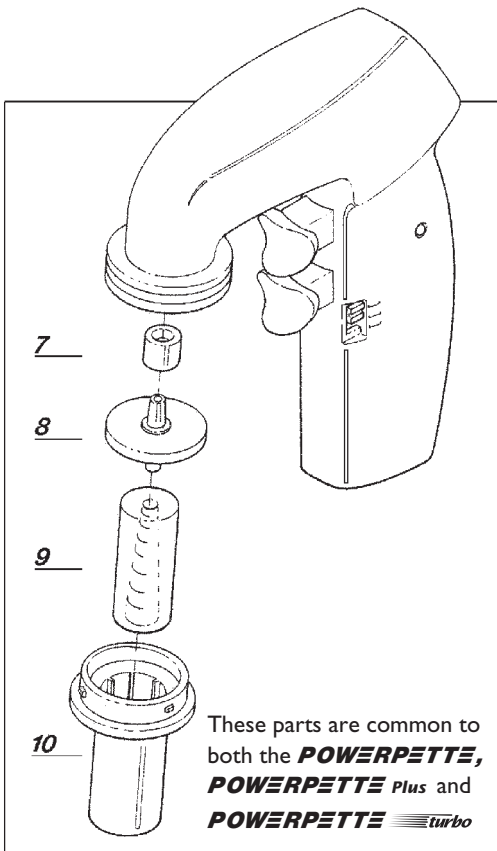Pipetting

To aspirate

- 1 Carefully hold the pipette and insert it securely into the polypropylene nose cone and silicone holder.
- 2 Select mode of operation – High, Low, or ‘G’.
- 3 Depress upper trigger (1) (The harder the pipetting triggers are pressed the faster the aspiration.)

To dispense

- 1 In BLOW-OUT mode, (mode switch on high or low) depress the lower trigger (2) (The harder the pipetting triggers are pressed, the faster the dispense).
- 2 If ‘G’ gravity mode is chosen, the upper aspirate trigger automatically operates at low speed while the lower trigger opens the pipette to atmosphere, allowing liquid to run out.

WARNING!

Please ensure that no liquid enters the instrument. The **POWERPETTE** / **POWERPETTE Plus** / **POWERPETTE turbo** should never be used without the filter (8). The filter also prevents liquid from entering the instrument. Where liquid has entered please clean and replace filter.

Trouble Shooting

TROUBLE	POSSIBLE CAUSE	REMEDY
Suction capacity slow	Filter blocked, dirty or damaged	Clean or replace membrane filter
	Batteries low	Recharge
Pipette drips	Pipette not securely positioned in silicone pipette holder	Re-attach
	Pipette holder or pipette damaged	Replace

The **POWERPETTE** / **POWERPETTE Plus** / **POWERPETTE turbo** are virtually service-free. If the recommendations above do not resolve the problem, please contact your authorized dealer.

Note

When the meniscus level has been reached liquid should not drip out of the pipette. If the pipette does drip, check that the adaptor assembly and filter are correctly assembled and the pipette is properly located into the silicone pipette holder.

Return for Repair

In the unlikely event of repair, or where the damage to the unit necessitates return to your distributor or the manufacturer, please decontaminate the unit and complete the required decontamination certificate where appropriate. Please also give details of the fault.

Voltages

UK	220V-240V 50/60Hz 3-pin
Europe	220V-240V 50/60Hz 2-pin
USA	100V-120V 50/60Hz 2-pin
Japan	100V-120V 50/60Hz 2-pin
India	220V-240V 50/60Hz 3-pin
Australia	220V-240V 50/60Hz 2-pin

Ordering Data

POWERPETTE complete with charger, bench stand and wall bracket with two spare membrane filters.

	UK	Europe	USA	Japan	India	Aust./NZ
	Cat. No.	Cat. No.	Cat. No.	Cat. No.	Cat. No.	Cat. No.
Grey	266-007	266-008	266-009	266-016	266-106	266-118
Yellow	266-069	266-070	266-071	266-072	266-107	266-119
Pink	266-073	266-074	266-075	266-076	266-108	266-120
Mauve	266-077	266-078	266-079	266-080	266-109	266-121
Blue/green	266-081	266-082	266-083	266-084	266-110	266-122
Mint green	266-085	266-086	266-087	266-088	266-111	266-123
Box of 5 spare 0.45 µm filters	266-003				
Box of 5 spare 0.20 µm filters	266-089				
0.45 µm sterile filter (1 off)	266-006				
Silicone Pipette holder	266-004				
Polypropylene nosecone	Please specify colour at time of order				

POWERPETTE Plus complete with charger, bench stand and two spare membrane filters.

	UK	Europe	USA	Japan	India	Aust./NZ
	Cat. No.	Cat. No.	Cat. No.	Cat. No.	Cat. No.	Cat. No.
Grey	266-000	266-001	266-002	266-015	266-100	266-112
Yellow	266-049	266-050	266-051	266-052	266-101	266-113
Pink	266-053	266-054	266-055	266-056	266-102	266-114
Mauve	266-057	266-058	266-059	266-060	266-103	266-115
Blue/green	266-061	266-062	266-063	266-064	266-104	266-116
Mint green	266-065	266-066	266-068	266-068	266-105	266-117
Box of 5 spare 0.45µm filters.....	266-003					
Box of 5 spare 0.20µm filters.....	266-089					
0.45µm sterile filter (1 off)	266-006					
Silicone Pipette holder	266-004					
Polypropylene nosecone	Please specify colour at time of order					

POWERPETTE turbo complete with charger, bench stand and two spare membrane filters.

	UK	Europe	USA	Japan	India	Aust./NZ
	Cat. No.	Cat. No.	Cat. No.	Cat. No.	Cat. No.	Cat. No.
Powerpette						
Turbo	266-130	266-131	266-132	266-133	266-154	266-160
Box of 5 spare 0.45µm filters.....	266-003					
Box of 5 spare 0.20µm filters.....	266-089					
0.45µm sterile filter (1 off)	266-006					
Silicone Pipette holder	266-004					
Polypropylene nosecone	Please specify colour at time of order					

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