

## Increase your productivity with ali-Q



**Faster aliquoting with less fatigue**

**Highly accurate and precise aliquots without the need to “eyeball” meniscus**



## Serological pipetting like you're used to...aliquoting like you've never been able to before!

ali-Q works just like any pipet controller but also stands like all Ovation® pipettes. However, unlike any other pipet controller, with ali-Q you can multi-dispense accurately and precisely with each push of the aliquoting button. You set the volume of each aliquot with the aliquoting dial, up to 5 mL.

### What sets Ovation ali-Q apart from every other standard pipet controller is its revolutionary aliquoting technology.

Simply turn the aliquot volume set dial to the volume you wish to repetitively dispense.

Then, aspirate as much volume as you will need to dispense the full number of aliquots. (Ex: You need to dispense 10 x 1 mL aliquots, then aspirate more than 10 mL)

Finally, press and hold the purple aliquot button to dispense each aliquot.

Eliminate time-consuming and inconsistent meniscus eyeballing! Let the ali-Q's state of the art intelligent measuring system do all the work to save you time and effort and increase your productivity!

#### VistaLab Catalog # 2000-0005

- Sensitive touch-control aspirate and dispense buttons
- Separate (purple textured) aliquoting button
- Aliquoting volume setting dial (purple)
- 0.45 µm replaceable membrane filter (1 extra included)
- Calibration certificate included (calibrated to Wobble-not serological pipet)
- Works with any size and brand of serological pipet
- Re-chargeable lithium ion battery (USB charger included)
- Wall mount included
- One-year warranty



For more info and to watch a video:  
<https://vistalab.com/ali-q-pipet-controller>

#### ali-Q features and benefits:

- Highly accurate and precise aliquoting without the need to "eyeball" the meniscus
- Push one button to aliquot
- Aliquot any volume from 0.5 to 5.0 mL with exceptional repeatability
- Compatible with all brands and sizes of serological pipets; optimal performance when using Wobble-not pipets
- Intelligent measuring system = real-time corrections for varying environmental conditions - No adjustment necessary!
- Accuracy and precision maintained at any dispense angle for more comfortable body positions, especially beneficial in the hood
- Ergonomic shape design for optimal grip comfort
- Self-standing, like all Ovation pipettes!