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## **Instruction Manual**

*Please read before use!*

# pipetus®

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Battery operated pipetting aid

P 990.70 01 / 1005

Hirschmann Laborgeräte  
GmbH & Co. KG  
Postfach 11 53 - Hauptstr. 7 - 15  
D- 74246 Eberstadt  
Germany  
Tel. +49(0)71 34/511-0  
Fax +49(0)71 34/511-90  
[www.hirschmann-laborgeraete.de](http://www.hirschmann-laborgeraete.de)  
[info@hirschmann-laborgeraete.de](mailto:info@hirschmann-laborgeraete.de)

- 2 -

## Inhalt

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## Contents

	Page
<b>1. Safety instructions</b>	5
<b>2. Components</b>	9
<b>3. Functional principle</b>	10
<b>4. Start-up</b>	12
4.1 Scope of delivery	12
4.2 Charging	13
4.3 Mounting the wall bracket	16
<b>5. Prohibited operations</b>	17
<b>6. Restricted operations</b>	18
<b>7. Pipetting</b>	19
7.1 Place the pipette	19
7.2 Fill the pipette	20
7.3 Empty the pipette	21
7.4 Operating speed adjustment	21

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<b>8. Spare parts replacement</b>	23
8.1 Replacement of pipetholder complete	23
8.2 Replacement of filter	23
8.3 Replacement of safety valve	24
8.4 Safety valve - Function check and troubleshooting	25
<b>9. Maintenance, cleaning and sterilisation</b>	27
9.1 Maintenance	27
9.2 Sterilisation	27
9.3 Cleaning	28
<b>10. Test of tightness</b>	29
<b>11. Repair at the manufacturer</b>	29
<b>12. Trouble shooting</b>	33
<b>13 Technical data</b>	35
<b>14. Accessories - spare parts</b>	37
<b>15. Clearance certificate form</b>	40

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*We are delighted that you have chosen our pipetting aid. The unit is a modern, extremely accurate pipetting aid which will facilitate your work considerably.*

*The pipetting aid is designed to the latest technical standards and is operationally safe. However, residual dangers can be associated with use of the unit, so it is imperative that the instruction manual be read carefully and all instructions therein heeded.*

*All the same, the instruction manual cannot illustrate every safety problem which may arise. Each user bears individual responsibility for adhering to safety and health regulations and determining and heeding limitations prior to using the unit.*

*Please observe the safety instructions prior to start-up in all cases!*

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## 1. Safety Instructions

- Observe all general safety practices and guidelines for laboratories, such as the use of protective clothing, safety goggles and gloves when handling hazardous chemicals.
- Never point the pipette in anyone's direction when in use!
- To avoid the risk of an electrical shock, never immerse either the unit or the recharger unit in water or any other liquid!
- Please ensure that the current of the power network matches with the voltage on the charger unit.

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- In the event of any damage of the cable or the recharger unit, the recharger unit can not be used.

- Never touch the plug of the recharger unit with wet hands!

- Observe the guidelines and procedures of the chemical manufacturers. (see also points 5 and 6)

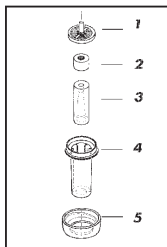
- The recharger unit may only be used in rooms. Protect from moisture!

- Non-rechargeable batteries should not be connected to the recharger unit!

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- A safety valve (2) in the pipetholder case (4) is located on top of the silicon adaptor (3) which prevents the liquid from getting into the handle, if the pipette is accidentally overfilled. If the safety valve has been activated, it must be removed by unscrewing the pipetholder and then cleaned and dried. Afterwards it must be inserted in the prescribed direction (see point 8). Wear safety gloves and goggles! (Danger of infection)

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**Attention!**

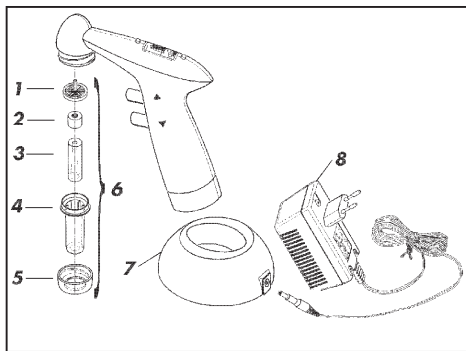
*If the unit is opened, the warranty is invalidated and any resultant claims void.*

- *The filter supplied contains a hydrophobic 0,2 µm membrane for air filtration in either direction.*
- *Only have the unit professionally serviced.*
- *Never use undue force!*

**Attention!**

*Only ever use the original recharger unit, accessories and replacement parts to avoid damaging the unit.*

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
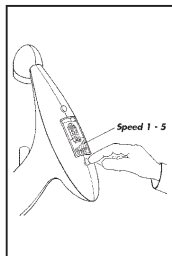
### 3. Functional principle

- The unit is a rechargeable pipetting controller for use with all standard pipettes from 0,1 ml to 200 ml.
- The pump is activated by pressing the In- ( $\Delta$ ) or respectively the Out- ( $\nabla$ ) button. The speed of filling (In) or dispensing (Out) is regulated by the intensity of pressure on the buttons.
- The pipetting speed is set by switching the 5 speed-stages and therefore an individual adaptation to the pipetted volume is possible.

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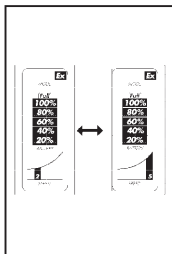
- The selected speed stage is indicated at the bottom of the LC display.

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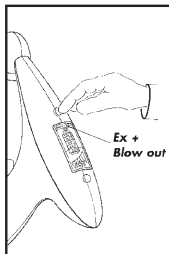- By switching from Ex to Blow out or vice versa, the unit can accommodate both to deliver and blow-out pipettes. In the Ex position the pump remains switched off while atmospheric pressure empties the pipette. In this way, the working time of the batteries is doubled.

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## 4. Start-up

### 4.1 Scope of delivery

The following components should be contained in the original carton:

- Pipetting aid
- 1 charging stand
- 1 power supply unit
- 1 wall bracket with 2 screws (chipboard-screws 3,0 x 30 mm) and 2 dowels (S5 \* 25 mm)
- 1 hydrophobic membrane filter 0.2 µm (yellow)
- 1 instruction manual

Please check that all parts have been delivered and are undamaged!

- The charger should not be used again if the cable or charger are damaged in any way!

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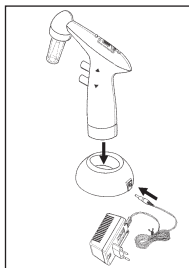
## 4.2 Charging

The unit is supplied with charged batteries.

However, the battery charge drops over time, and batteries should be charged prior to initial start-up.

- Either through inductive charging technology in the charging stand. Insert the plug of the charger unit into the socket of the charging stand.
- Or through direct charging of the unit. Insert the plug of the charger unit directly into the socket of the unit.
- Charging takes about 8 hours.

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### Guide:

The warming up of the unit during the charging operation is harmless.

### Attention!

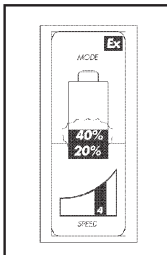
The unit may only be charged with the original recharger unit. The network voltage must match with the voltage indicated on the recharger unit (maximum deviation  $\pm 10\%$ ).

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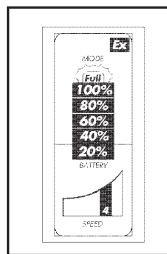
- One after the other segment of the battery capacity indication starts flashing during the charging procedure according to the respective charging status. Provided e.g. that the batteries are almost empty, the segment 20 % will be flashing.


- If the batteries are charged completely the display „Full“ will be flashing.

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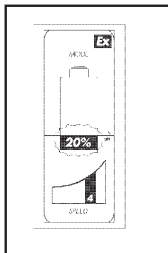- During pipetting one after the other segment of the display will disappear.
- When falling short of 10 % of the battery capacity, an acoustic signal will be heard after the pipetting operation.
- This indicates, that the unit must be charged again as soon as possible.

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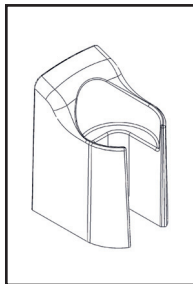
### 4.3 Mounting the wall holder

- Secure the wall holder to the wall in a suitable position with the two screws supplied (use the enclosed dowels if necessary). Drill two holes vertically positioned at a distance of 22 mm to each other for this purpose (drill size 5). Depth approx. 25 mm.
- Screw in the screws until a space remains between the wall and the screw heads.
- Suspend the pipette aid from the screws - Ensure that the unit rests lie flat to the wall.

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## 5. Prohibited operations

### **The unit can not be used with:**

- liquids of which vapours attack the materials silicone, PTFE, Novodur, PVDF, NBR, PCPA, PP and PA6.
- easily inflammable liquids (flame point below 0°C, for instance, ether and acetone)
- usage in explosive areas (Ex-area Zone 1)

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## 6. Restricted operations

- *Working and charging temperatures between +10°C and +40°C.*
- *Usage only in areas in which a dangerous, explosive atmosphere occurs only seldom and then only short-termed (Ex-area zone 2).*
- *Load only in Ex-secure areas.*
- *Do not keep unsupervised in an explosive area.*
- *liquids which emit copious volumes of gas, as vapours are deposited inside the pipetting aid and can corrode the unit.*

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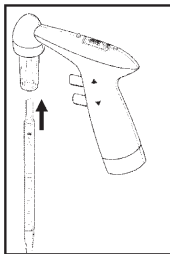
## 7. Pipetting

### 7.1 Place the pipette

- Place the pipette into the holder to ensure a secure and airproof fit!

#### **Attention!**

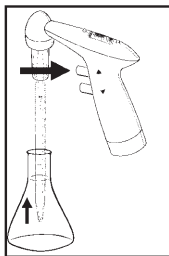
Danger of glass breakage, especially with thin pipettes.

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
## 7.2 Fill the pipette

- For filling the pipette, press the In-button ( $\Delta$ ).
- Select one of the 5 speed stages according to the volume of the pipette. Via the Min/Max button you can either chose minimum or maximum motor output (for larger volumina we recommend maximum output).

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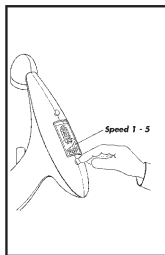### **Attention!**

No liquid should get into the unit as this can adversely affect the operating functions.

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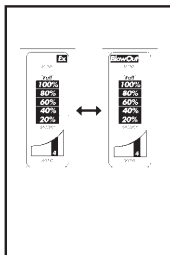
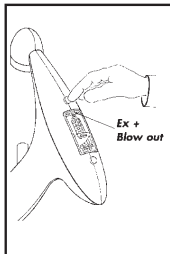
### 7.3 Empty the pipette

You can choose between two pipetting modes illustrated in the upper part of the display.


- Use of blow-out pipettes: If Blow out pipettes are used, the key has to be set so that „Blow out“ appears in the display yellow.
- ▶ Pipetting mode Blow out => pipette is blown out by the pump.
- Use of delivery pipettes: If To Deliver pipettes are used, the key has to be set so that „Ex“ appears in the display.
- ▶ Pipetting mode Ex => pipette is emptied by atmospheric pressure (without the pump).

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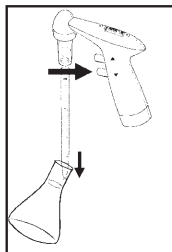=> To empty the pipette, press the Out button (▽), the intensity of finger pressure on the button regulates the working speed.

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#### 7.4 Operating speed adjustment

##### **Setting the suctioning speed:**

Press the In(△)button and simultaneously select the desired suctioning speed with the Speed button. By again pressing the In button, suctioning is done with the selected speed.

##### **Setting the blow-out speed:**

Press the Out(▽)button and simultaneously set the Speed button to the desired blow-out speed. The selected speed is then indicated through flashing.

The speed for the last mode selected (suctioning or blow-out) is always altered if the Speed button is pressed without simultaneously pressing the In or Out button.

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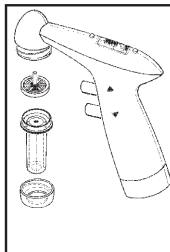
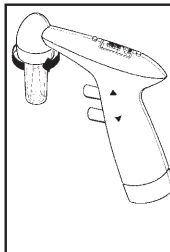
## 8. Spare parts replacement

### 8.1 Replacement of pipetholder complete

- Screw off the nut.
- Replace the pipetholder by a new one.

### 8.2 Replacement of filter

- Screw off the nut.  
Remove the old filter and place the new one in the pipetholder case so that the pivot shows upwards. Screw the pipetholder onto the unit, by turning the nut clockwise

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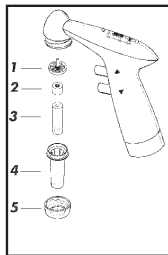
### 8.3 Replacement of safety valve

- Screw off the nut (5) counter-clockwise.

Take out the filter (1) and place it aside.

After that, remove the old safety valve (2) and place the new one in the pipetholder case (4), so that the black sealing ring shows upwards. Place filter (1)

on the safety valve with the pivot facing up. Screw the pipetholder onto the unit, by turning the nut clockwise.

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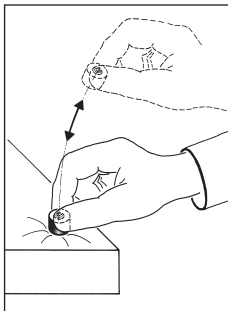
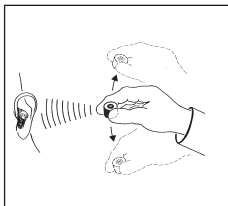
### 8.4 Function check and troubleshooting

#### IMPORTANT HINT FOR ASSEMBLING

Black O-Ring of safety valve must be visible upside above silicone adapter!

Function check of the safety valve. The membrane of the valve must create a "rattling" sound, while shaking the valve.

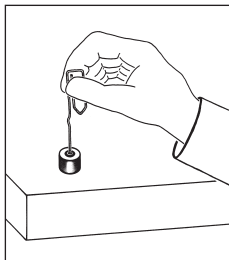
A sticking valve might become flexible again, by gently knocking it e.g. on the table edge. However in the event that the valve is glued, it must be soaked in a cleaning solution, rinsed with water and dried afterwards. Do not open the valve.

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
*In case the function of the valve is not restored yet, try to make it work again, by gently pushing the membrane with a thin and blunt object.*

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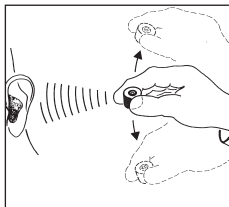## 9. Maintenance, Sterilisation and Cleaning

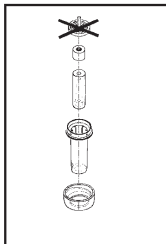
### 9.1 Maintenance

Under normal conditions of use, the unit is maintenance-free.

In the event, nevertheless, that they do become dirty or the suction becomes impaired, the safety valve and membrane filter should be replaced and the unit recharged.

### 9.2. Sterilisation

- The pipetholder, except the membrane filter, can be sterilised at 121°C.

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### 9.3 Cleaning

The safety valve (2) must be cleaned and dried before autoclaving. While reassembling the unit make sure that the side of the safety valve (2) with the black sealing ring shows towards the filter (1).

The pipetting aid handles can be cleaned with a damp cloth. Do not spray the unit, as liquid can penetrate to the interior.

#### **Attention!**

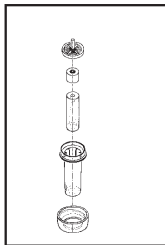
When cleaning and exchanging the safety valve and membrane filter, protective gloves and goggles must be worn (danger of infection)!

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## 10. Test of tightness

Once the required volume has been aspirated into the pipette, the liquid has to be kept in it unless the buttons are pressed. If the pipette however is dripping, disassemble the pipette holder and reassemble it again carefully.

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## 11. Repair at the manufacturer

Please observe the following points should it prove necessary to return the unit for repair (despite having handled it with all due care):

- Before sending the unit back, the user must ensure that it is carefully cleaned and decontaminated. The healthwise harmlessness must be confirmed, send the confirmation together with the unit to the manufacturer or the dealer. (A copy of the clearance certificate can be found on page 40)

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- Units will only be repaired, when they have been cleaned and decontaminated by the user.
- For safety reasons dirty units will be returned to the sender without being repaired.
- Please describe the malfunction exactly.
- Please return the unit complete with the power supply unit and charging stand in all cases, as this will facilitate checking of all components.
- Please quote the pipetted liquid.
- In the event of the unit being tampered with or being used in any improper way, (e.g. the application of undue force) any claims under the warranty are void.

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## 12. Trouble shooting

Symptom	Possible cause	Solution
<i>Liquid column in the pipette does not hold (pipette drops or runs out)</i>	<ul style="list-style-type: none"><li>- Safety valve built-in incorrectly</li><li>- Pipetholder is untight</li></ul>	<ul style="list-style-type: none"><li>- Insert the safety valve so, that the side with the black O-ring points to the filter (see point 8).</li><li>- Test the correct assembly sequence for the pipetholder. If required, test or replace individual components.</li><li>- Silicone adapters can become worn or cracked. Change if necessary</li><li>- Filters can be blocked. Change if necessary</li></ul>
<i>The pump will not blow out the pipette</i>	<ul style="list-style-type: none"><li>- Incorrect pipetting mode</li></ul>	<ul style="list-style-type: none"><li>- Set the Ex/Blow out key to the Blow out mode. The pipette will be blown out!</li></ul>
<i>The pump does not function neither filling nor blowing out</i>	<ul style="list-style-type: none"><li>- Batteries flat</li><li>- Safety valve incorrectly positioned</li></ul>	<ul style="list-style-type: none"><li>- Recharge the batteries with the recharger unit.</li><li>- Place the safety valve so that the black sealing ring shows towards the filter.</li></ul>

- 34 -

*Overfilled pipette*

*- Safety valve is activated immediately and prevents liquid from getting into the the unit*

*- Remove safety valve, clean and re-assemble. If required, replace the safety valve. During assembly take note of point 8.3.*

*Suction too slow or too weak*

*- Selected stage is too low*  
*- No tight, secure connection of the pipetholder*  
*- Filter is full*  
*- Silicone adaptor is damaged*

*- Improve the selection to a greater stage e.g. 3 -> 5.*  
*- Screw the pipetholder more tightly onto the handle.*  
*- Exchange the filter.*  
*- Exchange the silicone adaptor.*

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#### Batteries:

- 2 NiMH batteries
- Operation time for permanent use:
  - Position Ex 50 h
  - Position Blow out 25 h
- time for charging: 8 h

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*12 Monate Garantie*

CE:

*This mark is the confirmation that the unit conforms  
to the established requirements of the EU guidelines  
and has been tested according to these criteria.*

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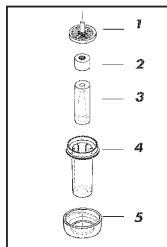
## 14. Accessories - spare parts

The spare parts for your pipetting aid listed below can be obtained from your specialised laboratory retailer:

Silicone adapter (3)	990 32 01
Safety valve (2)	990 32 02
Pipetholder akku complete	990 38 00
Pipetholder case (4)	990 38 03
Nut grey (5)	990 38 02
Round filter set (consisting of 5 replacement filters)	990 37 04
O-ring for cone	P990 32 05
Wall bracket (incl. dowels and screws)	990 56 00
Recharger unit complete with inductive charging stand, 230 V/50 Hz	990 75 03
Recharger unit complete with inductive charging stand, 240 V/50 Hz (UK version)	990 76 03
Recharger unit complete with inductive charging stand, 120 V/60 Hz	990 85 03

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..... (Serial number)

Medium used: .....

Blood, bodily fluids, patholog. samples:

**Yes**

**No**

**Please define in case of YES**

\_\_\_\_\_

Other hazardous biologic substances:

\_\_\_\_\_

Biologically-degradable materials which can pose a risk:

\_\_\_\_\_

Chemicals containing substances harmful to health:

\_\_\_\_\_

Radioactive materials:

\_\_\_\_\_

Other hazardous and toxic substances:

\_\_\_\_\_

### We hereby declare that

- the unit has been carefully cleaned and decontaminated prior to shipment

- no risks resulting from contamination exist

- we are aware that liability claims can be made against us in the event of damage being caused by contaminated units.

### Address of sender:

Name: \_\_\_\_\_

Date: \_\_\_\_\_

Signature: \_\_\_\_\_