Transferpette®

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Operating Manual

Please read carefully before use,
and follow all operating and safety
instructions!

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Zur Reparatur einsenden

Entsorgung

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| Seite | Contents | Page |
|--------------|---|-------------|
| 4 | Safety Instructions | 4 |
| 4 | Purpose | 4 |
| 5 | Limitations | 5 |
| 5 | Operating Elements | 5 |
| 6 | Handling | 6 |
| 6 | - Use proper pipette tips | 6 |
| 6 | - Volume Setting | 6 |
| 6 | - Hand posture while pipetting | 6 |
| 7 | - Ejecting the tip | 7 |
| 8 | Pipetting | 8 |
| 8 | - Aspirating the sample | 8 |
| 9 | - Dispensing the sample | 9 |
| 10 | Maintenance | 10 |
| 10 | - Lubricating the plunger | 10 |
| 11 | Adjustment | 11 |
| 11 | - Fixed-volume models | 11 |
| 12 | - Variable-volume models | 12 |
| 14 | Checking the volume | 14 |
| 15 | Calculation of accuracy/coefficient of variation | 15 |
| 16 | Cleaning | 16 |
| 18 | Autoclaving | 18 |
| 21 | Troubleshooting | 22 |
| 23 | Specifications / Spares | 23 |
| 26 | Accessories | 26 |
| 28 | Return to repair | 28 |
| 28 | Disposal | 28 |
| 29 | Warranty | 29 |

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Read carefully before use!

Safety Instructions

1. Observe general hazard prevention instructions and safety regulations; e.g. wear protective clothing, eye protection and gloves.
2. Observe the reagent manufacturer's information.
3. Read this Operating Manual carefully.
4. Always use the instrument in such a way that neither the user nor any other persons are endangered.
5. Only use manufacturer's original parts.
6. Operate only with tip attached.
7. Use original PLASTIBRAND® pipette tips, or make sure that the tips to be used are compatible.

Purpose

Air-displacement pipette for pipetting aqueous solutions of medium density and viscosity.

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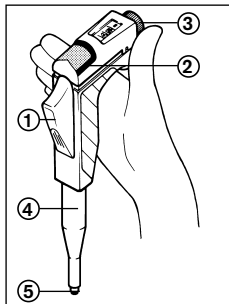
Operating Limitations

The user has to ensure the compatibility of the instrument with the intended application.

This instrument cannot be used:

- for liquids incompatible with polypropylene,
- for liquids of a very high vapour pressure,
- in Reverse Mode (ISO 8655/2).

Viscous and wetting liquids may impair the accuracy of the volume. The same applies to liquids of a temperature deviating from the ambient temperature by more than $\pm 5\text{ }^{\circ}\text{C}$ (permissible operating range: $15 - 40\text{ }^{\circ}\text{C}/59 - 104\text{ }^{\circ}\text{F}$).

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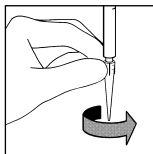
Handling

- Use proper pipette tips
Pipette performance can only be guaranteed if high-quality pipette tips are used. We consequently recommend the use of PLASTIBRAND® pipette tips.
- pipette tips are disposables


Fitting the tip

- in regular intervals, check the pipetting system - consisting of Transferpette® and pipette tip - for tightness.
- turn them lightly in the direction of the arrow

Note:

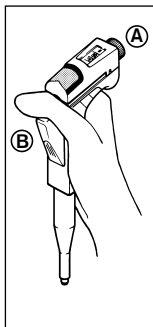If the tip is turned in the other direction, the screw connection of the shaft tip may be loosened, with air leaks resulting. The shaft tip must be tightened at the latest when a red ring becomes visible on it.

(A) Volume setting

(B) Hand posture while pipetting

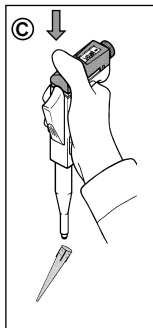
- Put your thumb across the pipetting key
- unlike with conventional pipettes.

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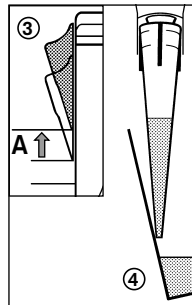
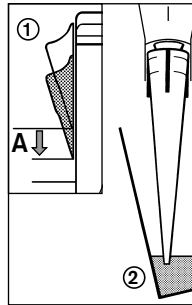
Pipetting

Aspirating the sample

- Rinse the tip once with the sample liquid.

1. Press the lateral pipetting key to **the first stop (A)**.
2. Immerse the pipette tip 2 - 3 millimeters into the sample.

Note:

Leave the tip immersed in the liquid for about another second to avoid the intake of air.

3. Let the pipetting key slide back **slowly**.
4. Lightly wipe the tip against the wall of the vessel.

Note:

Don't lay the instrument horizontal when the tip is full, or liquid may penetrate inside the instrument.

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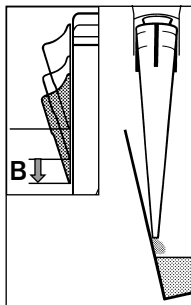
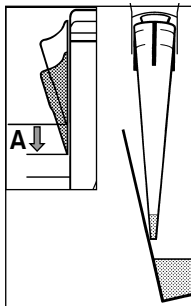
Dispensing the sample

Place the pipette tip against the wall of the vessel.

- Press the pipetting key down to the stop (A) and hold it down.

Note:

For serum and liquids of high viscosity or low surface tension, observe adequate waiting time to improve accuracy.

- The blow-out stroke empties the tip completely: Press the pipetting key down to the stop (B).

Note:

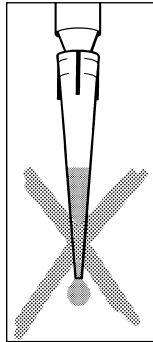
When working with serum and liquids of high viscosity, low surface tension or at low volumes: Rinse the tip with the reagent contained in the vessel to improve the accuracy.

- Wipe the pipette tip against the wall of the vessel.
- Let the pipetting key slide back.

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Maintenance

In regular intervals:

1. Check firm seat of pipette shaft cone, tighten if necessary.
2. Examine piston and seal for contamination (see page 16 "Cleaning").
3. Check volume.

Lubricating the plunger

Very thinly lubricate the plunger:

- after cleaning,
- if motion is uneven,
- after autoclaving.

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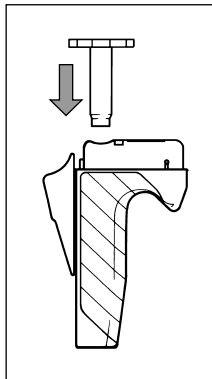
Adjustment

The instrument is permanently adjusted for aqueous solutions. It can be adjusted for solutions of other density and viscosity.

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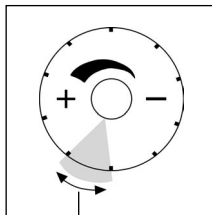
1/250 vom Nennvolumen

Note:

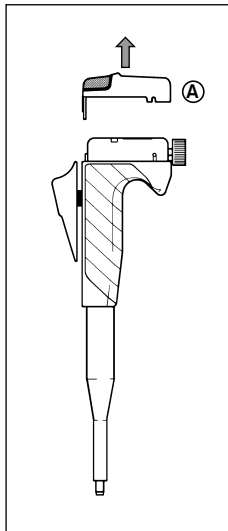
Maximum permissible adjustment is 3 full turns.

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
Variable-volume models


1. Check the volume, determine actual value (see page 14, "Checking the volume").
2. Release lateral closures by turning into position .
3. Pull off ejector cap (A).

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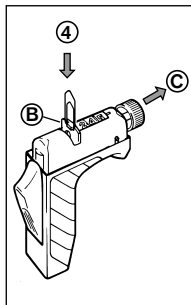4. Use a paper clip to extract the red distance plate (B) from the top.
5. Pull out volume setting knob (C) to the stop in the direction of its axis (approx. 2 mm).

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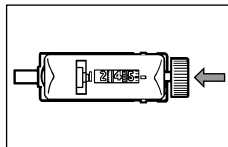6. Set to the previously measured actual value. Push back volume setting knob to the stop (approx. 2 mm), without turning it.
7. Replace the red distance plate.
8. Mount ejector cap (A).
9. Secure ejector cap by turning the lateral closures into position \ominus .

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Checking the volume

The gravimetric volume test of the pipette requires the following steps:
(Testing procedures are described e.g., in ISO 8655 Part 6).

Adjust volume to nominal capacity.

Conditioning the pipette:

To condition the pipette before testing, mount a new pipette tip and pipette five times with testing liquid (H_2O dist.). Discard the pipette tip.

Carrying out the test (observe instructions by the balance manufacturer):

- Mount a new pipette tip and pre-rinse once with testing liquid.
- Take up testing liquid and pipette into the weighing vessel.
- Weigh the pipetted quantity with an analytical balance.
- Calculate the volume, taking the temperature into account.
- At least 10 pipettings and weighings in three volume ranges (100%, 50%, 10%) are recommended for statistical analysis.

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Variationskoeffizient

$$\text{VK\%} = \frac{100 s}{\bar{V}}$$

Calculation of accuracy (A%) and coefficient of variation (CV%):

Accuracy (A%) and coefficient of variation (CV%) are calculated according to the equations used in statistical quality control.

Calculations (for the nominal volume)

$$\text{Mean value } \bar{x} = \frac{\sum x_i}{n} \quad \begin{array}{l} x_i = \text{results of weighings} \\ n = \text{number of weighings} \end{array}$$

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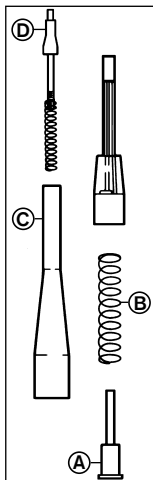
Cleaning

Note:

The Transferpette® requires regular maintenance, depending on frequency of use and pipetted media.

After cleaning, always check volume (see page 14).

Pipettes up to 200 µl

(Grey and yellow colour code)

1. Eject the tip.
2. Unscrew pipette shaft (C) from the handle.
3. Remove plunger (A) and plunger spring (B) from the handle.
4. Unscrew pipette shaft cone (D) from the pipette shaft (C).
5. Clean parts shown on drawing with soap solution or isopropyl alcohol. Afterwards rinse with distilled water.
6. Dry the parts (max. 120 °C/248 °F).
7. Lubricate plunger very thinly ("Maintenance", page 10).
8. Reassemble parts after they have cooled down.

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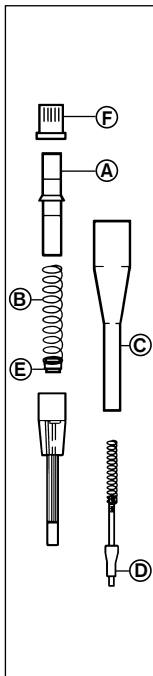
Pipettes of 200 µl upwards

1. Eject the tip.
2. Unscrew pipette shaft (C) from handle.
3. Unscrew retention bushing (F) from pipette shaft (C).
4. Remove plunger (A) and plunger spring with seal (B), (E).

Note:

Do not disassemble parts (B) and (E) any further. Take care not to deform the seal; clean it only with a soft lint-free cloth.

5. Unscrew pipette shaft cone (D) from the pipette shaft (C).
6. Clean parts shown on drawing with soap solution or isopropyl alcohol. Afterwards rinse with distilled water.
7. Dry the parts (max. 120 °C/248 °F).
8. Lubricate plunger ("Maintenance", page 10).
9. Reassemble parts after they have cooled down.

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– continued –

Note:

Do not overtighten the retention bushing (F). Tighten only by hand.

10. After assembling, operate the pipetting lever five times to centre the seal.

Autoclaving

For sterile work or for decontamination, the pipette parts shown here withstand steam sterilization at 121 °C/250 °F (2 bar absolute) for 20 min. exposure time (t_e), according to DIN.

Note:

To protect against contamination, filter tips can also be used.

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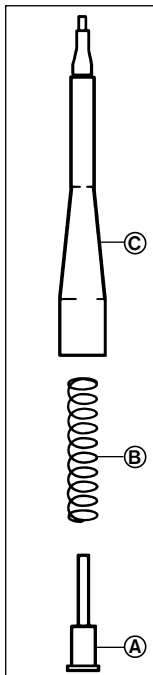
**Attention:**

It is user's responsibility to ensure effective autoclaving.

After each autoclaving:

- Check volume (see page 14).

Note:

Only autoclave parts that have been cleaned.

Pipettes up to 200 µl**(Grey and yellow colour code)**

1. Eject the tip.
2. Cautiously unscrew pipette shaft (C) from handle.
3. Remove plunger (A) and spring (B) from the handle.

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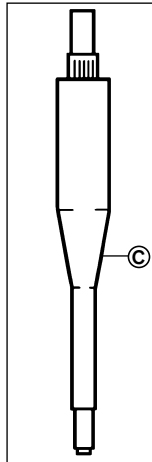
Pipettes up to 200 μl

- continued -

4. Sterilize the complete shaft, plunger and plunger spring in the autoclave.
5. Allow the parts to cool completely and assemble in the reverse order.

Pipettes of 200 μl upwards

1. Eject the tip.
2. Cautiously unscrew pipette shaft (C) from handle.
3. Autoclave the complete shaft (C) in the autoclave without having it disassembled.
4. Allow to dry and to cool completely.
5. Screw pipette shaft into handle.

Note:

The handle is not autoclavable. For disinfection you may use commercially available disinfecting solutions as long as they do not attack PP.

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Troubleshooting

| Trouble | Possible Causes | Action to be taken |
|---|--|--|
| <ul style="list-style-type: none">- Tip dripping (instrument leaks)- Aspirated volume too low- Delivered volume too low | <ul style="list-style-type: none">- Unsuitable tip- Tip not seated tightly- Pipette shaft cone loose - Seal contaminated- Seal damaged- Non-aqueous solutions pipetted- Piston damaged | <ul style="list-style-type: none">- Only use high-quality tips- Press tip on firmly- Tighten clockwise see "Handling", page 6see "Cleaning", page 16see "Specifications/Spares", page 25see "Adjustment", page 11see "Spares", page 25 |
| <ul style="list-style-type: none">- Slow suction or none | <ul style="list-style-type: none">- Shaft clogged | <ul style="list-style-type: none">- Clean shaft see "Cleaning", page 16 |
| <ul style="list-style-type: none">- Volume too high | <ul style="list-style-type: none">- Pipetting key pressed too far (into the blow-out position) before sample take up. | <ul style="list-style-type: none">- Operate properlysee "Pipetting", page 8 |

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Specifications/Spares

(Spare parts can be fitted by the user)

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* Error limits relative to the nominal capacity (maximum volume) indicated on the instrument, obtained at equal temperature (20 °C) of instrument, ambience and dist H₂O, and with smooth, jerk-free operation. The absolute error margins (µl) are not exceeded at any partial volume.


** Corresponding to the color of the tips.

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Following tolerances are typically (manufacturer's experience) obtained by new instruments under optimized testing conditions (qualified operators and standardized ambience conditions):

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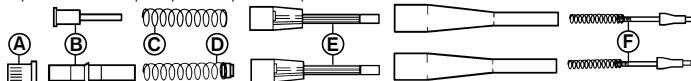*** Pipette shaft complete with parts (A) to (F)

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Accessories

| | |
|---|------------------|
| Rack 3 instruments, | Cat. No. 7032 03 |
| Rack 2 x 3 instruments, | Cat. No. 7032 08 |
| Wall support for 3 instr., | Cat. No. 7032 10 |
| Silicon oil V 140 000 mm ² /s, | Cat. No. 7032 05 |
| Spring (C) | Cat. No. 9884 |
| Retention bushing (A) | Cat. No. 6530 92 |

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Note:

ISO 9000 and GLP-guidelines require regular examinations of your volumetric instruments. We recommend to check the volume every 3 - 12 months. The interval depends on the specific requirements on the instrument. For instruments frequently used or in use with aggressive media, the interval should be shorter.

The detailed testing instruction can be downloaded on www.brand.de or can be ordered under Cat. No. 9976 19 (english) in printed form.

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Return for Repair

Attention:

Transporting of hazardous materials without a permit is a violation of federal law.

BrandTech Scientific, Inc. will not accept instruments that are not appropriately cleaned and decontaminated.

- Therefore, contact BrandTech Scientific, Inc. and obtain return authorisation **before** sending your instrument for service.

Disposal

For the disposal of instruments and tips, please observe the disposal regulations valid in your area.

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Warranty

As provided by law, any and all warranties are null and void if the product has been misused, modified or repaired by unauthorized personnel, if the defects are caused by negligence (instruction manual, maintenance) or by normal wear and tear. Use only original manufacturer's accessory/spare parts.

For conditions and extent of warranty refer to our General Conditions of Sale (www.brandtech.com).

Subject to technical modification without notice. Errors excepted.

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© Spitze abwerfen
BrandTech Scientific, Inc. wird keine Geräte annehmen, die nicht ordnungsgemäß gereinigt und dekontaminiert wurden.